X17C SAFETY CHECKLIST					BROOKHAVEN NATIONAL LABORATORY NATIONAL SYNCHROTRON LIGHT SOURCE
Rev: B	Effective: 5/24/07		Page 1 of 1		Number: LS-SCL-0040
Reviewed by:	Thomas McDonald	Reviewed by	y :		Approved by: J.Aloi

Original signatures on file

The only official copy of this file is the one on-line in the NSLS Quality Assurance website. Before using a printed copy, verify that it is the most current version by checking the document effective date on the NSLS QA website.

	1.	Valid Padlock Index and log, no locks open
	2.	Cable penetration shield upstream of double plug doors in place (top of wall - see photo)
	3.	Cement bricks in place along entire bottom length of the two X17 front end sawtooth plug doors
		(see photo)
	4.	Water cooled backstop in place in downstream end of X17C hutch (see photo).
		Bremsstrahlung Shield BS6 in place and banded (see photo from item #4).
		Note: BS3, BS4, BS5 are part of X17B checklists
	6.	Bremsstrahlung Shield BS1 (in front of MS Monochromator) in place and banded (see photo)
		Inside Transport Hutch (view items through window if hutch is interlocked)
	7.	Bremsstrahlung Shield BS2 (downstream end of hutch) in place and banded (see photo)
		Inside Transport Hutch (view items through window if hutch is interlocked)
	8.	Threshold shield in place under door on X16 side of hutch
		Threshold shield in place under door on X18 side of hutch
		X17B transport pipe lead shielded and strapped.
		Upstream X17B transport pipe wall penetration lead shields in place (see photo)
		Downstream X17B transport pipe wall penetration lead shields in place (see photo from item #4
Ω	DC() / Reamline Staff: