## Detection of CYP1A1 in Formalin-Fixed, Paraffin-Embedded Rat Tissue

## **Reagent and Antibody Information**

1X Wash Buffer
3% Hydrogen Peroxide
1% BSA Diluent
1X Citrate Buffer
DAB Chromagen
Hematoxylin

Staining Kit: Vectastain Elite ABC Kit (Rabbit IgG)
Vector Laboratories, Inc.
Burlingame, CA 94010
www.vectorlabs.com
1-800-227-6666
Catalog # PK-6101

**Note**: This kit contains all reagents necessary to make the blocking reagent, secondary antibody and label complex.

Avidin / Biotin Blocking Kit Vector Laboratories, Inc. Burlingame, CA 94010 www.vectorlabs.com 1-800-227-6666 Catalog # SP-2001

Primary Antibody: Rabbit Anti-Rat CYP1A1 Polyclonal Antibody

Chemicon International, Inc. Temecula, CA 92590 www.chemicon.com 1-800-437-7500

Catalog # AB1247

Negative Control Serum: Normal Rabbit Serum Jackson Immunoresearch Laboratories, Inc.

West Grove, PA 19390

www.jacksonimmuno.com

1-800-367-5296

Catalog # 011-000-001

## **Staining Procedure**

Positive Control Tissue: Pancreas (islets of Langerhans) and TCDD-treated liver Stain Localization: Cytoplasmic (centrilobular pattern in the liver)

1. Deparaffinize and hydrate slides through the following solutions:

Solution	Repetitions	Time
Xylene	2 times	5 minutes
100% Ethanol	2 times	3 minutes
95% Ethanol	2 times	3 minutes
1X Wash Buffer	2 times	5 minutes

- 2. Quench endogenous peroxidase by placing the slides in 3% hydrogen peroxide for 15 minutes.
- 3. Rinse the slides in 2 changes of 1X Wash Buffer for 5 minutes each.

	Vectastain Rabbit Elite Staining Kit Exp Date New Kit: yes / no
1.	Apply the block from the Rabbit Elite Kit, and incubate for 20 minutes at room temperature.
	DO NOT RINSE THE SLIDES. CONTINUE TO AVIDIN-BIOTIN BLOCK.
5.	Avidin / Biotin Blocking Kit  Lot # Exp. Date New Kit: yes / no  Apply avidin block for 15 minutes at room temperature.  Quick rinse in 1X Wash Buffer.  Apply biotin block for 15 minutes at room temperature.
	DO NOT RINSE SLIDES WITH BUFFER BEFORE ADDING PRIMARY ANTIBODY.

6. Apply primary antibody at a 1:100 dilution, and incubate for 1 hour at room temperature.

Lot #\_\_\_\_\_\_ Exp Date \_\_\_\_\_\_

For negative control slides, dilute the protein concentration of the normal rabbit serum to match that of the primary antibody. Make a 1:100 dilution from this normalized serum, and apply to the slides. Incubate for 1 hour at room temperature.

Lot #\_\_\_\_\_ Date Reconstituted\_\_\_\_\_

7. Rinse the slides in 2 changes of 1X Wash Buffer for 5 minutes each.

ONLY WIPE EXCESS BUFFER.

- 8. Apply the secondary antibody from Rabbit Elite Kit, and incubate for 30 minutes at room temperature.
- 9. Rinse the slides in 2 changes of 1X Wash Buffer for 5 minutes each.

- 10. Apply the label complex from the Rabbit Elite Kit, and incubate for 30 minutes at room temperature. (Prepare at least 30 minutes prior to use.)
- 11. Rinse the slides in 2 changes of 1X Wash Buffer for 5 minutes each.

12. Apply the DAI	3 chromagen, and incubate in	the dark for 6 minutes	at room temperature.
(Add 1 drop of	DAB per ml of substrate)		
Lot #	Exp. Date	New Kit:	yes / no

- 13. Rinse the slides in tap water 3 minutes.
- 14. Counterstain with Harris Hematoxylin for 20 seconds.
- 15. Rinse the slides in tap water until water is clear.
- 16. Gently agitate slides in 1X Wash Buffer until they turn blue.
- 17. Dehydrate through the following solutions:

Solution	Repetitions	Time
95% Ethanol	1 time	3 minutes
100% Ethanol	3 times	3 minutes
Xylene	2 times	5 minutes

18. Coverslip

Updated 10/04/04