X11A SAFETY CHECKLIST				]	BROOKHAVEN NATIONAL LABORATORY NATIONAL SYNCHROTRON LIGHT SOURCE
Rev: C	Effective: 5/24/2010		Page 1 of	1	Number: LS-SCL-0024
Prepared by: Philip Marino		Reviewed by: P. Marino			Approved by: B. Chmiel

Original signatures on file.

The only official copy of this file is the one on-line in the NSLS Quality Assurance website. Before using a printed copy, verify that it is the most current version by checking the document effective date on the NSLS QA website.

□ 1.	Valid Padlock Index and log, no padlocks open.
	Exclusion Zones (EZ #1) and (EZ #2) in place and secure covering Vacuum Bellows #1 wrapped in lead, Acrylead on Viewport, Water cooled Beryllium window #1 wrapped in lead, 8" Conflate flange covered with lead sheet.
□ 3.	Bremsstrahlung shield (BS #1) in place and banded as per photo
	Lead scatter shielding as per photo
□ 5.	Exclusion Zones (EZ #3) and (EZ #4) in place and secured.
	Viewport #1 and #2 on top of monochromator covered with lead glass and lead sheet. (Note: Bremsstrahlung shields BS # 2 and BS #3 under lead, not visible),
□ 7.	Lead shielding above the monochromator tank is in place as per photo.
□ 8.	Lead shielded mono venting box in place.
□ 9.	Two layers of acrylead on top Viewport, one layer on side Viewport, and 5 way cross covered in lead (near the photon shutter).
OPC	O / Beamline Staff: Date:
	User System Checks
	Beam-line isolation valve 1A is in the open position. Vacuum gauge on front of monochromator reads 28" or higher.
Beam	line Rep.: Date: