X17B1 SAFETY CHECKLIST					BROOKHAVEN NATIONAL LABORATORY NATIONAL SYNCHROTRON LIGHT SOURCE
Rev: C	Effective: 5/24/2010		Page 1 of 1		Number: LS-SCL-0037
Prepared by: Philip Marino		Reviewed by: P. Marino			Approved by: B. Chmiel

Original signatures on file.

The only official copy of this file is the one on-line in the NSLS Quality Assurance website. Before using a printed copy, verify that it is the most current version by checking the document effective date on the NSLS QA website.

- $\hfill\square$ 1. Valid Padlock Index and log, no locks open
- \Box 2. Cable penetration shield upstream of double plug doors in place as per photo.
- □ 3. Cement bricks in place along entire bottom length of the two X17 front end sawtooth plug doors (see photo)
- □ 4. X17B1 hutch exterior (right of door) lead-covered KF blank flange over vacuum port when not in use.
- □ 4A. X17B1 hutch exterior (X17C side) lead-covered KF blank flange over vacuum port when not in use.
- □ 5. X17B1 hutch interior water cooled backstop/ Bremsstrahlung shield #3 down stream end of hutch (see photo)
 - Banded and in the beam path, Kirk key removed and placed in X17B1 Backstop SRU (outside hutch next to safety shutter control panel) for X17B1 operation

OR

- Banded and out of the beam path for X17B2/B3 operation
- □ 6. Bremsstrahlung Shield #1 (in front of MS Monochromator) in place and banded (see photo **Inside Transport Hutch** (view items through window if hutch is interlocked))
- □ 7. Bremsstrahlung Shield #2 (downstream end of hutch) in place and banded (see photo) **Inside Transport Hutch** (view items through window if hutch is interlocked)
- □ 8. X17B1 hutch threshold shielding bolted in place under door going into hutch.

OPCO / Beamline Staff: Date:	
------------------------------	--