

GENERAL INSTRUCTION FOR DOING WIPE TESTS:

Enclosed are envelopes containing number wipes.

- (1) Fill in the blanks on the envelopes with the information requested (i.e. location and ECD source number).
- (2) After the information is written on the envelopes, put on the glove.
- (3) Remove one wipe, thoroughly rub around the exhaust port of the Electron Capture Detector (ECD) and put the wipe back into the envelope. Repeat this procedure with the other wipes for each detector.
- (4) Discard the glove and return the envelopes as soon as possible.